

P1 Experiment
Detection

EXP

11 am

Subj. #: 307

Date: 6-2-97

Subject:

Subj. type: Focal Lesion

Amp. Settings (make note of any changes)

	Gain	HPF	LPF
Ch. 1-28	.10	.01	100
Ch. 29	.10	.01	100
Ch. 30	.10	.01	100
Ch. 31	.10	.01	100

External
filter

50

Run Calibration

On DIG:

```
cd \dig
digitize sf.hdr <filename>.log <filename>.raw 1
```

On STIM:

```
cd \experiments\p1
cal
c 990
s [on DIG]
s [on STIM]
```

After approximately 30 cal pulses,

```
p [on DIG]
q [on STIM]
```

Cap placement

mid/ro

Head circumference: _____ cm

Nasion-to-inion: 34 cm

Channel Assignments: (make note of any irregularities)

1: FPz	9: FC2	17: CP1	25: PO7
2: FC (plug-in) RUC	10: FC6	18: CP2	26: PO3
3: F3	11: T7	19: CP6	+27: POz
4: Fz	12: C3	20: P7	+28: PO4
5: F4	+13: C4	21: P3	29: PO8
6: FC (plug-in) LC	14: Cz +	22: Pz	30: O1
+7: FC5	15: T8	23: P4	31: Oz
8: FC1	16: CP5	+24: P8	32: O2

Comments: